


Pasteur Pipette Adapter and Flow Controller

Pasteur Pipette Adapter and Flow Controller

Two types of Pasteur pipette adapters are offered.

Pasteur Pipette Adapter and Flow Controller, Male Luer to Fluted Nut - NA0636


A. Loosen the fluted nut

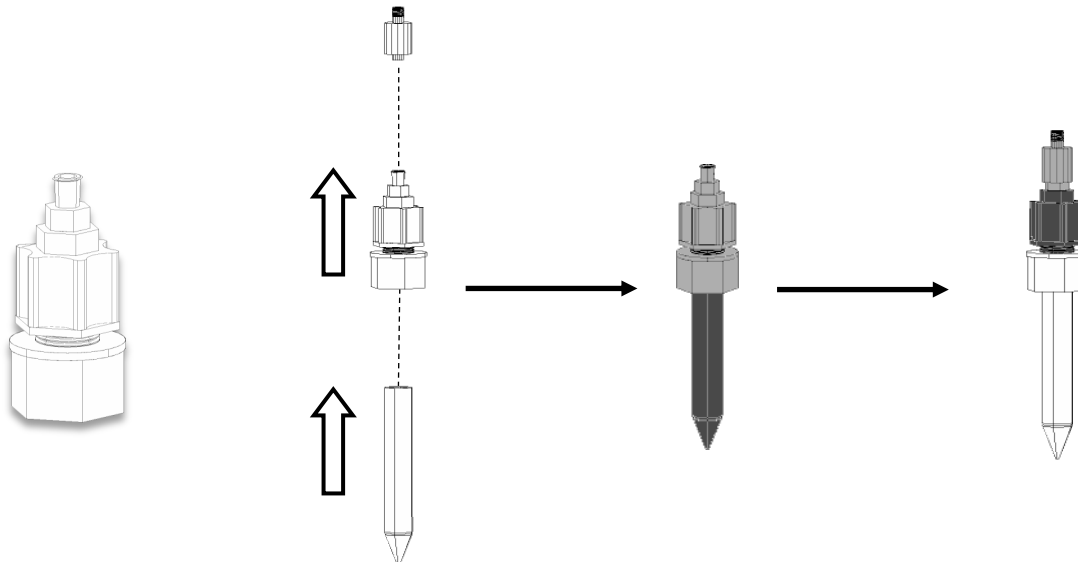

B. Push pipette into the fluted nut and tighten.

C. Screw the adapter onto the luer fitting like you would a needle.

Pasteur Pipette Adapter and Flow Controller, Male Luer to 1/8 Hose Barb - NA0637

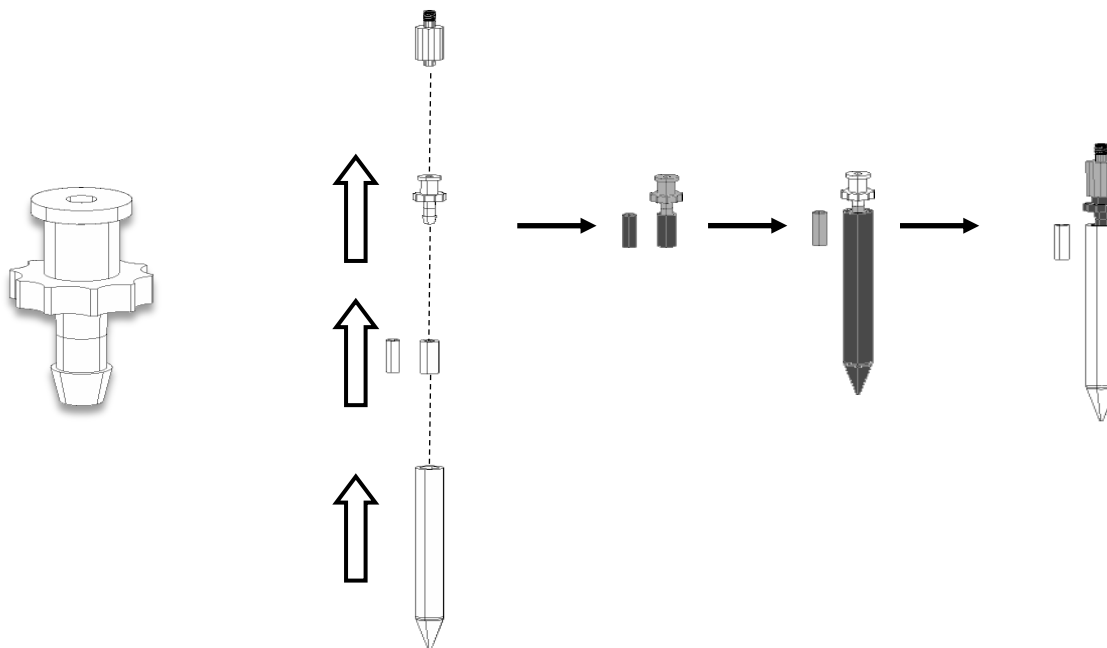

A. Choose the tube that fits your pipette.

B. Put the tube on the end of the adapter.

C. Push the pipette onto the tube on the adapter.

D. Screw the adapter onto the luer fitting like you would a needle.